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Alarm Code Table

Code No.	Cause	Remedy
01	Attempt to start turning the rotor with the door opened	(1)Close the door completely.
02	In a state of standby for low degree vacuum for 30 minutes or more	(1)Open the door and check for complete closing. (2)If the evaporator is moistened,press DEFROST
03	The vacuum switch is not actuated upon lapse of 5 minutes or more after the vacuum pump is turned ON.	Ditto
04	The vacuum switch is turned OFF at 5,000rpm or more.	(1)Eliminate the cause of deterioration of vacuum.
05	Temperature above the limit.	(1)If the vacuum has been teriorated,eliminate its cause.
06	No rotation signal from the rotor.(Operation without a rotor)	(1)If no rotor is set in the centrifuge,set one in it. (2)The adapter, if defective,should be replaced with a normal one.
07	Operation with the rotor speed presetat a level over the limit.	(1)Lower the presetting.
08	Abnormal rotation signal from the rotor	(1)Clean the overspeed adapter of the rotor. (2)Replace the overspeed adapter with a new one
09	Unbalance of the rotor is detected.	(1)Eliminate the cause of the unbalance
10	Above the speed limit of rotor.	(1)Check to see if the preset rotation speed is not above the limit.
11	Circuit breaker of the lubricant feed pump (oil pump) turned off.	(1)Contact a service technician for instruction on repair.
12	Rotor started after power failure (however,the alarm code "12" is displayed when the speed at power recovery is lower less than 1,000rpm than the speed setting)	(1)Centrifugation condition are variable.Make an alteration if necessary.
13	Deceleration in response to a signal from an accessory unit.	(1)The cause lies in an accessory unit.Check the accessory units.
20	Abnormal rotation signal from motor.	(1)If no rotor set in the centrifuge,set one in it.
81	Change the memoried operation condition.	(1)After press the reset switch reset the operation condition.(including the memory of step mode or etc.)